

FocalCheck™ Fluorescence Microscope Test Slides

F36909 FocalCheck™ fluorescence microscope test slide #1

F36913 FocalCheck™ fluorescence microscope test slide #2

F36914 FocalCheck™ fluorescence microscope test slide #3

Quick Facts

Storage upon receipt:

- Room temperature

applications. The microspheres are mounted in optical cement (refractive index of ~1.52) for maximal stability and are intended for use on all fluorescence microscope systems.

- **FocalCheck™ fluorescence microscope test slide #1** is ideal for routine checking and calibration of confocal and widefield fluorescence microscopes.
- **FocalCheck™ fluorescence microscope test slide #2** provides a robust, reproducible method of evaluating the performance of spectral imaging systems, as well as the ability to discriminate closely overlapping spectra. The slide consists of 6 µm-diameter microspheres labeled with a series of spectrally overlapping dyes.
- **FocalCheck™ fluorescence microscope test slide #3** is useful for basic evaluation of filter performance and as a general practice slide for fluorescence microscopy and digital imaging.

Introduction

FocalCheck™ fluorescence microscope test slides #1, #2, and #3 are specifically designed for calibrating fluorescence microscope systems and evaluating system and filter performance. The slides each contain 10 sample areas coated with proprietary fluorescent microspheres designed specifically for microscopy

Table 1. Contents of FocalCheck™ fluorescence microscope test slide #1.

	1	2	3	4	5
Row A	15 µm	6 µm	4 µm	1 µm	0.5 µm
	Green/orange/far red thin rings	Blue throughout/green ring	Blue/green/red/far red	Blue/green/red/far red	Blue/green/red/far red
Row B (Intensity Series)	6 µm	6 µm	6 µm	6 µm	6 µm
	Green	Green	Green	Green	Green
	100% intensity	33% intensity	10% intensity	3% intensity	0.667% intensity

Table 2. Contents of FocalCheck™ fluorescence microscope test slide #2.

	1	2	3	4	5
Row A	6 µm	6 µm	6 µm	6 µm	Mixture of all four double-stained beads
	Green	Orange	Red	Far red	
	Bead 1 503/511 nm Bead 2 511/524 nm	Bead 1 541/555 nm Bead 2 545/565 nm	Bead 1 578/605 nm Bead 2 589/613 nm	Bead 1 657/676 nm Bead 2 671/692 nm	
Row B	6 µm	6 µm	6 µm	6 µm	Mixture of all eight reference beads
	Green	Orange	Red	Far red	
	Ring 503/511 nm Core 511/524 nm	Ring 541/555 nm Core 545/565 nm	Ring 589/613 nm Core 578/605 nm	Ring 671/692 nm Core 657/676 nm	

Table 3. Contents of FocalCheck™ fluorescence microscope test slide #3.

	1	2	3	4	5
Row A	2.5 μm	2.5 μm	2.5 μm	2.5 μm	2.5 μm
	Blue	Green	Orange	Red	Deep red
	350/440 nm	505/515 nm	540/560 nm	580/605 nm	633/660 nm
	Bright	Bright	Bright	Bright	Bright
	100% intensity	100% intensity	100% intensity	100% intensity	100% intensity
Row B	2.5 μm	2.5 μm	2.5 μm	2.5 μm	2.5 μm
	Blue	Green	Orange	Red	Deep red
	350/440 nm	505/515 nm	540/560 nm	580/605 nm	633/660 nm
	Dim	Dim	Dim	Dim	Dim
	1% intensity	1% intensity	1% intensity	1% intensity	1% intensity

Storage and Handling

The FocalCheck™ fluorescence microscope test slides should be stored at room temperature. Avoid prolonged exposure to light. When stored correctly, the slides should be stable for at least 12 months.

Protocol

The basic procedure for using the FocalCheck™ fluorescence microscope test slides is to image the beads using standard techniques. Below are several general applications.

Color Registration

For multicolor experiments, particularly on widefield microscopes, switching between different filter configurations can result in misregistration between the various color channels. Using the multicolor beads on FocalCheck™ fluorescence microscope test slide #1 (sample positions A1–5), the extent of the misregistration can be digitally corrected independently of the sample (Figure 1).

Chromatic Distortions

Modern microscopes use highly color-corrected (achromatic) optical components. However, it is possible that certain optical components may introduce wavelength-dependent image shifts that can affect the correct registration of multicolor labels. Such aberrations are critical to avoid in colocalization experiments. FocalCheck™ fluorescence microscope slide #1 is ideal for ensuring that the imaging system is free from chromatic distortions.

X- and Y-Axis Calibration

For distance and other morphometric measurements, microscope imaging should be calibrated for each objective and magnification setting using a stage micrometer. The larger 15 μm beads on FocalCheck™ fluorescence microscope test slide #1, sample positions A1 and A2, provide ideal targets for assessing the calibration of the microscope in three dimensions. When measured at the equator, the beads are nominally 15 μm in diameter; this should correspond with the measured X and Y dimensions by imaging.

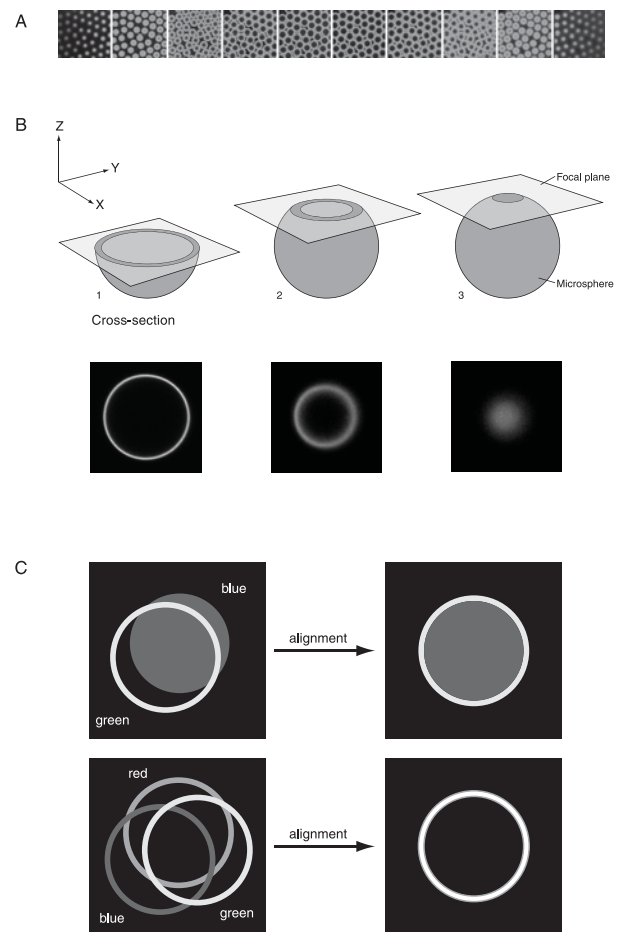


Figure 1. Confocal laser-scanning microscope optical cross-sectioning and alignment with FocalCheck™ microspheres. A) Serial optical sectioning from top to bottom along the z-axis of ring-stained microspheres reveals a continuous pattern of disc-to-ring-to-disc images. B) The diameter of the fluorescent ring (or disc) seen is dependent on the depth of the optical focal plane. C) In the confocal laser-scanning microscope, separate light paths exist for UV and visible wavelengths. Also, emitted fluorescence is detected by different photomultipliers. Proper optical alignment may be obtained with either of two types of FocalCheck™ microspheres. For example, the microspheres with a green ring stain that are blue-fluorescent throughout the bead allow UV/visible wavelength alignment in three dimensions upon aligning the green ring with the blue disc. Focal alignment is also possible simultaneously in three colors by aligning the red, blue, and green rings of the FocalCheck™ microspheres containing fluorescent red, blue and green ring stains.

Z-Axis Calibration

Positions A1 and A2 on FocalCheck™ fluorescence microscope test slide #1 can also be used to assess the performance of the motorized focus control of an imaging system. To do this, acquire Z-series stacks from bottom-to-top and top-to-bottom. The refractive index of the slide mounting media is close to 1.515, which minimizes spherical aberration when using an oil immersion lens. When viewed in orthogonal projections, the bead images should appear relatively circular. Any distortions could indicate problems in the Z-stepping of the microscope’s focus system or a refractive index mismatch.


Lamp Stability

Most standard mercury lamps are rated for 200–300 hours of use. While many can last longer, there is often little warning that the lamp is failing. Bead samples on any of the FocalCheck™ fluorescence microscope test slides can be used to monitor the stability of the excitation source on the microscope. Periodic imaging of the same sample, using the same acquisition parameters, can indicate changes in lamp output by plotting the intensity of the beads over time.

Intensity Calibration

FocalCheck™ fluorescence microscope test slide #1 also contains a series of intensity beads. By acquiring a series of images, you can create a relative intensity curve against which samples can be evaluated (Figure 2). The intensity of the beads can be used to generate curves for different instruments, thus providing a basis for assessing the relative fluorescence intensity of a sample in a system-independent manner.

FocalCheck™ fluorescence microscope test slide #1 can also be used to examine the linearity and dynamic range of the detection system on the microscope under various settings. The intensity range of the beads was chosen to cover the approximate dynamic range of a typical 12-bit CCD camera (4096 gray levels) when imaged at a single exposure setting.

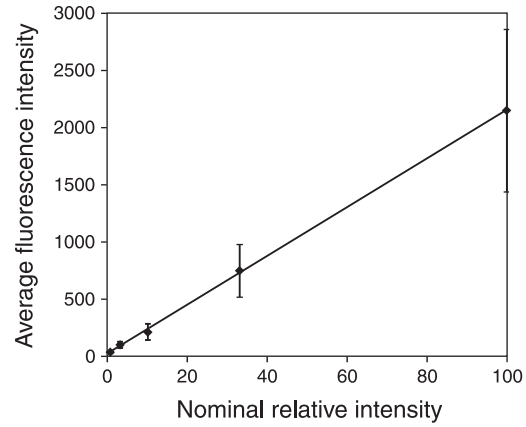

Nominal relative intensity	1.00%	5.00%	10.00%	40.00%	100%
Actual intensity	18.81	88.70	201.91	738.01	2138.48
Determined relative intensity	0.88%	4.15%	9.44%	34.51%	100%

Figure 2. Average fluorescence intensity of the intensity beads series on FocalCheck™ microscope test slide #1. The intensity bead series was imaged on a Nikon Eclipse 800 microscope using a 100 watt Hg-arc lamp and a 60x/1.4NA objective. The images were acquired using a Hamamatsu Orca-ER 12-bit CCD camera. The average fluorescence intensity and standard deviation of the bead targets was determined and plotted against nominal relative intensity as determined by flow cytometry. In this particular configuration, all five intensities were imaged using a single exposure setting. The measured average intensities shown in the table give relative intensities that correspond closely to the nominal values measured by flow cytometry.

Spectral Scanning

FocalCheck™ fluorescence microscope test slide #2 can be used on spectral imaging instruments to help ensure that the wavelength detection system is working correctly. This slide can also be used to perform many of the general tests described above.

Product List *Current prices may be obtained from our Web site or from our Customer Service Department.*

Cat #	Product Name	Unit Size
F36909	FocalCheck™ fluorescence microscope test slide #1 *for alignment, intensity, and calibration*	each
F36913	FocalCheck™ fluorescence microscope test slide #2 *for spectral imaging systems*	each
F36914	FocalCheck™ fluorescence microscope test slide #3 *5 colors, high and low intensities*	each
P7220	PS-Speck™ Microscope Point Source Kit *blue, green, orange and deep red fluorescent beads*	1 kit

Contact Information

Further information on Molecular Probes products, including product bibliographies, is available from your local distributor or directly from Molecular Probes. Customers in Europe, Africa and the Middle East should contact our office in Paisley, United Kingdom. All others should contact our Technical Service Department in Eugene, Oregon.

Please visit our website — probes.invitrogen.com — for the most up-to-date information.

Molecular Probes, Inc.

29851 Willow Creek Road, Eugene, OR 97402
Phone: (541) 465-8300 • Fax: (541) 335-0504

Customer Service: 6:00 am to 4:30 pm (Pacific Time)

Phone: (541) 335-0338 • Fax: (541) 335-0305 • probesorder@invitrogen.com

Toll-Free Ordering for USA:

Order Phone: (800) 438-2209 • Order Fax: (800) 438-0228

Technical Service: 8:00 am to 4:00 pm (Pacific Time)

Phone: (541) 335-0353 • Toll-Free (800) 438-2209

Fax: (541) 335-0238 • probestech@invitrogen.com

Invitrogen European Headquarters

Invitrogen, Ltd.

3 Fountain Drive

Inchinnan Business Park

Paisley PA4 9RF, UK

Phone: +44 (0) 141 814 6100 • Fax: +44 (0) 141 814 6260

Email: euroinfo@invitrogen.com

Technical Services: eurotech@invitrogen.com

Molecular Probes products are high-quality reagents and materials intended for research purposes only. These products must be used by, or directly under the supervision of, a technically qualified individual experienced in handling potentially hazardous chemicals. Please read the Material Safety Data Sheet provided for each product; other regulatory considerations may apply.

Limited Use Label License

For research use only. Not intended for any animal or human therapeutic or diagnostic use. The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) to not transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of the above patents based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Molecular Probes, Inc., Business Development, 29851 Willow Creek Road, Eugene, OR 97402. Tel: (541) 465-8300. Fax: (541) 335-0504.

Several Molecular Probes products and product applications are covered by U.S. and foreign patents and patents pending. All names containing the designation ® are registered with the U.S. Patent and Trademark Office.

Copyright 2005, Molecular Probes, Inc. All rights reserved. This information is subject to change without notice.